

## Release notes – Clemex Software v9.0.314.25775

Release Date: 01-11-2021

### Bug Fixes

| Product                   | Description   | Notes |
|---------------------------|---|-------|
| Powder Disperser          | New recognize the new pressure gauge model. (940)   |       |
| Properties Administration | Clear Last Input now work properly. (923)   |       |
| CMT                       | Stage window is now in foreground focus when selected. (887)  |       |
| CIR                       | EN10247 report is now available. (909)  |       |
| CIR                       | DataBrowser : ASTM E45-18 E Results format, data browser do not show same particles for B Thin and B Heavy anymore. (907)         |       |
| CIR                       | ISO 4967 analysis, Export Raw data in txt and CSV is now working properly. (854)  |       |
| VisionPE/CIR/CMT          | Camera switching on Nikon MA 200 is now working properly. (845)   |       |
| CIR                       | Normal mode - Export Results to CSV is now working properly. (771)  |       |
| Vision                    | Tofra LED Driver Deadlock When Closing Vision is resolved. (939)  |       |
| Vision                    | Toolbox Instructions: Validation Issue - keyboard entry, Tab no longer exit field resolved. (881)                                 |       |
| Vision                    | Properties: "Field Count" is now displayed in correct format. (471)   |       |
| Vision                    | Toolbox - Binary Operations – operation parameters have now the same range limits using the keyboard and the up down arrow. (374) |       |
| Vision                    | Load images instruction behaviors are now fixed. (643)  |       |

### New Features

| Product     | Description                                      | Notes   |
|-------------|--|---------|
| Vision      | Martensitic Toolbox Instruction version 2.0      | UPGRADE |
| Vision      | Austenitic Toolbox Instruction v1.0              | NEW     |
| Vision      | Perlite-Ferrite Toolbox Instruction v1.0         | NEW     |
| App Package | "Martensitic" application package                | NEW     |
| App Package | "Austenitic" application package                 | NEW     |
| App Package | "Perlite-Ferrite" application package            | NEW     |
| App Package | "Aluminium/Magnesium Alloys" application package | NEW     |

**Known Issues/Bugs** (to be fixed in a later release)

| Product |  | Workaround  |
|---------|--|---|
| Vision  | Window size scaling issue when new instructions “Martensitic/Austenitic/Perlite-Ferrite” are used. (959)       | Set OS windows size scaling to: 100%                                  |
| Vision  | Object measure issue, ordering of measurement items may freeze the app at run time. (960)                      | None  |
| Vision  | “Martensitic/Austenitic/Perlite-Ferrite” instructions parameter values may not be set by keyboard entry. (964) | Use the value field Up/Down arrow control to change parameter values. |