

Release notes – Clemex Software v9.2.xxx.xxxx RTM

Release Date: ??

Software Bug Fixes

#	Product	Description	Notes
VPES-1148	User Management	AuditViewer3 must not allow to delete backed up audit trail databases	
VPES-1124	User Management	Audit Management must not be facultative	
VPES-1143	Captiva	Manual Multi-Layer Grab HDR processing must not include an Auto-Exposure as the user must have the ability to do this by himself	
VPES-1082	PSA	Auto-exposure issue with camera Clemex 2.1M USB2 (03-221)	
VPES-1104	Vision PE/Lite	Calibration properties in report, Tolerance issue	
VPES-1110	Vision PE/Lite	Special text annotation do not work properly in other language than English.	
VPES-1111	Vision PE/Lite	Auto Calibration German translation issue - Calibrate All Mags	
VPES-1118	Vision PE/Lite	Vision prevents automatic Windows screen lock	
VPES-365	Vision PE/Lite	Video Input Operation - Many Operation parameters, input shall not accept floating point or scientific notation due to space constraint.	
VPES-1097	Vision PE/Lite	Fix Character Set Issues in Tracing Code	
VPES-362	Vision PE/Lite	Autofocus per lens settings - multiple minor entries validation issues	
VPES-366	Vision PE/Lite	Video Input Operation - Build Mosaic, Max Mosaic Size parameter does not appear to be validated.	
VPES-368	Vision PE/Lite	Toolbox - Frame Operations - Insert Guard Frame, multiple validation issues	
VPES-867	Vision PE/Lite	Focus offset and focus interval are not updated automatically	
VPES-910	Vision PE/Lite	Incorrect ASTM E112 grain size menu	
VPES-1117	Vision PE/Lite	No response when using imported Plugins	
VPES-1129	Vision PE/Lite	Routine containing ClemexStudio plugin cannot be saved	
VPES-1154	Vision PE/Lite	Strange Hang when running Plugin instruction	Todo: paul
VPES-990	PSFilter	Toolbox 'Undo' operation doesn't work properly	
VPES-867	CMT	Focus offset and focus interval are not updated automatically	
VPES-1054	Clemex Software Setup	Implement Auto Deinstallation of Previous version of Clemex Soft from the install program	In progress: paul
VPES-1152	Clemex Software Setup	CMT v9.1 cannot be installed alone - A Firewall Plugin instruction is generated.	Todo: paul
VPES-1027	CIR	CIR Max width ISO4967 Issue	In progress: Sebastien

New Features

#	Product	Description	Notes
VPES-1108	User Management	New "Act as an Auditor" built-in General protected component	In progress: Sebastien
VPES-980	Vision PE / Lite / CMT / CIR / Captiva	Now Windows 11 Compatible	There is multiple issues found in RC1 that need to be fixed: Sebastien

Hardware Support

#	Issue	Brand	Device Type	Model

Known Issues/Bugs (to be fixed in a later release)

#	Product		Workaround
VPES-958	Vision PE/Lite	In some rare situation, disconnecting from Excel report engine may crash the application.	Reboot the PC and retry.
VPES-676	Vision PE/Lite	Errors when using 'X/Y projection' measures, value(s) of the X and/or Y projection may be wrong on the label.	none
VPES-1015, VPES-764	Vision PE/Lite	In some situation building mosaic can crash the application	Reboot the PC and retry.
VPES-990	Vision PE/Lite	Toolbox 'Undo' operation doesn't work properly	Change value back manually
VPES-982	CIR	Run workspace standard selection combo box is not always behaving properly under win11	none
VPES-870	CIR	map edition - cannot transfer to/from Beta for EN10247	none
VPES-436	CIR	Load Image Mode - It shall not be possible to edit heat and sample	none
VPES-439	CIR	Normal mode - switching camera also activates detection but buttons are off.	none
VPES-513	CIR	Open stage file issue - if a new and unsaved pattern is present, opening an existing pattern will merge the 2.	none
VPES-1163	CMT	Stage pattern window shortcut issues.	none