

Release notes – Clemex Vision PE/Lite v9.1.628.26016.LT9 RTM

Release Date: March 10th, 2022

Software Bug Fixes

#	Product	Description	Notes
VPES-960	Vision PE/Lite	Ordering of measurement items do not freeze the application anymore – (960)	
VPES-959	Vision PE/Lite	No more windows sizing issues caused by new Plugins UI (959)	
VPES-925	Vision PE/Lite	Square grid instruction is now working properly (925)	
VPES-965	Vision PE/Lite	Toolbox - IF System Condition, second condition is now set in routine (925)	
VPES-998	Vision PE/Lite	Chart condition is now working properly (998)	
VPES-964	Vision PE/Lite	Plugin Instructions parameters can now be set by keyboard entry (964)	
VPES-364	Vision PE/Lite	Auto Exposure, now accept upper limit by keyboard input. (364)	
VPES-361	Vision PE/Lite	Rectangular and Elliptical Field Template fields validation issues fixed	
VPES-974	ConfigSaver	ConfigSaver utility is now working normally (974)	
VPES-744	User Management	Workstation info is now available from audit trail (774)	
VPES-829	User Management	Changing user password now reset the countdown for expiration (829)	

New Features

#	Product	Description	Notes
VPES-992	Vision PE	Latest Micrometer certification updated (992)	

Hardware Support

#	Issue	Brand	Device Type	Model
VPES-977	Bug	Leica	Stereoscope	M205A – Is now fully compatible (977)
VPES-993	New	Leica	Stereoscope	M205C – Initial Integration (993)
VPES-994	New	Leica	Stereoscope	M165C – Initial Integration (994)

Known Issues/Bugs (to be fixed in a later release)

#	Product		Workaround
VPES-990	Vision PE/Lite	New plugin instructions cannot be undone with Toolbox Undo button	Modify the instruction parameters manually.